STEAM-THRU® SERIES CONNECTORS

Steam-Thru® Connectors allow a quick and easy sterile connection between stainless steel biopharmaceutical processing equipment and disposable bag and tube assemblies. The single-use design saves time and money by eliminating unnecessary cleaning procedures and reducing validation burden associated with reusable components.

SPECIFICATIONS

OPERATING CONDITIONS (Fluid Transfer)

STEAM POSITION

Temperature: 39°F to 104°F (4°C to 40°C) Pressure: Up to 30 psi, 2.1 bar (Steam-Thru) Up to 35 psi, 2.4 bar (Steam-Thru II)

FLOW POSITION

Temperature: 39°F to 104°F (4°C to 40°C) **Pressure:** Vacuum to 20 psi, 1.4 bar

STERILIZATION

 Full Connector Assembly
 Gamma: Up to 50 kGy irradiation
 Autoclave: Up to 265°F (129°C) for 60 minutes
 Steam-In-Place (SIP):
 Up to 266°F/130°C (up to ~24psi),
 60 minutes (Steam-Thru)

Up to 275°F/135°C (up to ~31psi), 60 minutes (Steam-Thru II)

TERMINATIONS

3/8" (9.5mm) to 1/2" (12.7mm) ID hose barb (Steam-Thru) 3/8" (9.5mm) to 1/2" (12.7mm) ID hose barb and 3/4" sanitary (Steam-Thru II)

MATERIALS

Connection: Polysulfone (amber tint) **O-rings:** Silicone (clear), platinum-cured **Removeable Sleeve:** Polycarbonate

NOTE: Steam pressures are estimated based upon information in Steam Tables found in literature



BENEFITS

Allows a true steam-through SIP process which eliminates "dead legs" and the need for laminar flow hoods

Allows sterile connection and disconnection and permits high media flow rate

Secures valve position, provides visual indicator of process stage $% \left({{{\bf{r}}_{\rm{s}}}} \right)$

Easily connects to process equipment

TYPICAL FLOW RATE:

Innovative three-port design

Steam-Thru II thumb latch

3/4" and 1-1/2" sanitary terminations

Patented valve design

Cv Value Range:

FEATURES

4.0 - 7.0 for Steam-Thru 3.0 - 9.0 for Steam-Thru II



Cv values represent the approximate expected flow rate in gallons per minute of water at room temperature for a 1 PSI pressure drop. The flow is generally constrained by the smallest diameter, which in some cases will be the termination diameter and not the Nominal Flow Path.

Scan code to visit webpage



cpcworldwide.com/STC

NOTE

Validation and Extractables data can be requested at cpcworldwide.com/STC

DID YOU KNOW

Steam-Thru connectors are perfect for any hybrid processing at your facility. If using a stainless-steel bioreactor and single-use systems, easily make a sterile connection between the two systems with a Steam-Thru connector on your single-use system and mounting it directly onto your bioreactor.



COLDER PRODUCTS COMPANY

+1 (651) 645-0091 +1 (651) 645-5404 EE: +1 (800) 519-7633

PHONE-

COLDER PRODUCTS COMPANY GMBH Germany PHONE: +49 6026 9973-0
 COLDER PRODUCTS COMPANY LIMITED

 ShangHai, China

 PHONE:
 +86 21 2411 2666

 TOLL FREE:
 +86 400 990 1978

STEAM-THRU SERIES DIMENSIONS


POLYSULFONE with polycarbonate sleeve



TERMINATION PART NO. A B G D e 3/4" X 3/4" SANITARY X 1/2" HB STC1700000 1.20" (30.5 mm) 5.09" (129.3 mm) 4.44" (112.8 mm) 2.00" (50.8 mm) 0.89" (22.6 mm) 3/4" X 3/4" SANITARY X 3/8" HB STC1700100 4.80" (121.9 mm) 4.15" (105.4 mm) 2.00" (50.8 mm) 0.60" (15.2 mm) 1.20" (30.5 mm) 3/4" X 1-1/2" SANITARY X 1/2" HB STC1700200 1.20" (30.5 mm) 5.09" (129.3 mm) 4.44" (112.8 mm) 2.00" (50.8 mm) 0.89" (22.6 mm) 3/4" X 1-1/2" SANITARY X 3/8" HB STC1700300 4.15" (105.4 mm) 1.20" (30.5 mm) 4.80" (121.9 mm) 2.00" (50.8 mm) 0.60" (15.2 mm)

STEAM-THRU CONFIGURATIONS

Steam-Thru Connection's patented three-port design allows steam to pass directly through the lower ports to "steam on" to stainless equipment. After the SIP cycle is completed, the connector's valve is actuated, creating a sterile flow path to single-use systems.

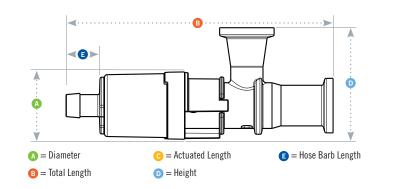

STEAM-THRU II SERIES DIMENSIONS

POLYSULFONE with polycarbonate sleeve



TERMINATION	PART NO.	A	B	C
3/4" X 3/4" SANITARY X 1/2" HB	STC2020000	1.42" (36.1 mm)	6.84" (173.7 mm)	5.93" (150.6 mm)
3/4" X 3/4" SANITARY X 3/8" HB	STC2020100	1.42" (36.1 mm)	6.76" (171.7 mm)	5.93" (150.6 mm)
3/4" X 1-1/2" SANITARY X 1/2" HB	STC2020200	1.42" (36.1 mm)	6.84" (173.7 mm)	5.93" (150.6 mm)
3/4" X 1-1/2" SANITARY X 3/8" HB	STC2020300	1.42" (36.1 mm)	6.76" (171.7 mm)	5.93" (150.6 mm)
3/4" X 3/4" SANITARY X 3/4" Sanitary	STC2020900	1.42" (36.1 mm)	6.60" (167.6 mm)	5.93" (150.6 mm)

All measurements are in inches (millimeters) unless otherwise noted. Tubing must meet stated inside and outside diameters.

2.40" (61.0 mm)

Ø

0.88" (22.4 mm)

0.80" (20.3 mm)

0.88" (22.4 mm)

0.80" (20.3 mm)

0.62" (15.7 mm)


0.62" (15.7 mm)

STEAM-THRU II CONFIGURATIONS

3/4" X 1-1/2" SANITARY X

3/4" SANITARY

Steam-Thru II Connections offer the flexibility of "steam on" and "steam off" functionality. The innovative design allows the valve to be returned to the steam position enabling a second SIP cycle following media transfer. The "steam off" disconnection of single-use systems minimizes cross-contamination risks associated with reusable components.

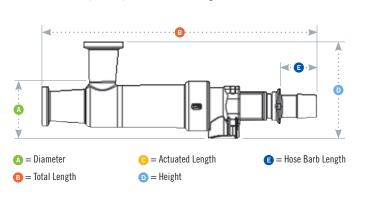

WARRANTY: All sales are subject to Colder Products Company's limited express warranty set forth in the CPC catalog. Contact your local distributor or CPC Customer Service for warranty provisions. Warning: Due to the wide variety of possible fluid media and operating conditions, unintended consequences may result from the use of this product, all of which are beyond the control of CPC. It is the user's responsibility to carefully determine and test for compatibility for use with their application. All such risks shall be assumed by the buyer.

STC2021000 1.42" (36.1 mm) 6.60" (167.6 mm) 5.93" (150.6 mm)

COPYRIGHT © 2021 BY COLDER PRODUCTS COMPANY. COLDER PRODUCTS COMPANY, COLDER PRODUCTS AND CPC ARE REGISTERED

TRADEMARKS WITH THE US PATENT & TRADEMARK OFFICE